

From Eye to Insight



## QUICK GUIDE

### Enersight software

How to update Emspira 3 digital microscope and Flexacam C3 microscope camera



1

Download the Firmware file for Emspira 3 or Flexacam C3 and copy it on a USB stick. They can be found at:

Emspira 3: <https://www.leica-microsystems.com/products/digital-microscopes/p/emspira-3/>

Flexacam C3: <https://www.leica-microsystems.com/de/produkte/mikroskop-kameras/p/flexacam-c3/>

2

Plug the USB stick into your Emspira 3 or Flexacam C3

3

Turn on the on-screen display (OSD) - if not done yet

4

Navigate to Settings > General

5

Click on "Update" button underneath "Current Version"

6

The system restarts automatically and it's ready to use!

